





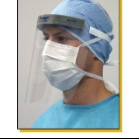


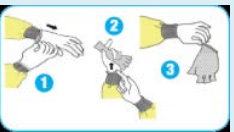


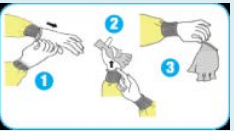






PUTTING ON (DONNING) Personal Protective Equipment

1	Wash hands	
2	Change into scrubs (if required)	
3	Put on booties (if required)	
4	Put on Gown or Lab coat	
5	Put on Head cover (if required)	
6	Put on face mask (if required) <ul style="list-style-type: none"> N-95 respirator for infectious (must be fit tested annually) Surgical mask for animals 	
7	Put on eye protection or face shield (if required)	
8	Put on gloves – use extended cuff length if available. If 2 nd pair of gloves is required, the extended 12” cuff is preferred for the top layer.	
While wearing PPE avoid touching your face or adjusting the PPE once in the infectious area.		

REMOVING (DOFFING) Personal Protective Equipment

You are removing PPE from the most contaminated to least.

1	Remove booties (if wearing)	
2	Remove gloves <ul style="list-style-type: none"> If wearing double gloves remove outer gloves only 	
3	Remove gown or lab coat	
4	Wash hands or remove 2 nd pair of gloves	 
5	Remove eye protection or face shield (if wearing)	
6	Remove face mask <ul style="list-style-type: none"> Handle straps only, do not touch the front portion 	
7	Remove head cover (if wearing)	
8	Wash hands	
9	Change into street clothes (if wearing scrubs) and wash hands	